

GLBIOCONTROL

R A P I D M I C R O B I A L D I A G N O S T I C

OPERATING MODE DENDRIDIAG® BF

Measuring the quantity of bacteria on surfaces by ATP-metry

- MANAGEMENT OF CLEANING & DISINFECTION -

CONTENT

GL BIOCONTROL	page 4
What is quantitative ATP-metry?	page 5
Applications	page 6
Equipment needed	page 7
Protocol	page 8
Interpretation of results	page 11
Troubleshooting	page 13
Controls	page 14
Contact	page 15

GL BIOCONTROL

GL BIOCONTROL specializes in environmental risk management and is an expert in sanitary engineering along with biological monitoring of water and surfaces.

Our clients are environment professionals: industrialists, industry operators, water treatment companies, laboratories, study design engineer and public authorities.

We have several skills including: risk assessments, research and development, analysis, product development and professional training. Through these areas of expertise, GL BIOCONTROL:

- **Develops** risk management tools (ATP-metry kits for total flora quantification in water, surface or air, concentration kit for virus analyses...).
- **Uses** methodologies and innovative tools to study the microbial world (qPCR, NGS, ATP-metry...).
- **Studies** ecosystems to anticipate and prevent public health risks, in particular linked to *Legionella* and *Pseudomonas* genus (risk assessment, microbiology diagnostics, ATP cartography...).
- **Advise**s water sector professionals on how to manage their facilities in order to reduce public health risks as well as improve the environmental footprint (water, treatment products and energy conservation).
- **Trains** environmental professionals on microbiological risk management and laboratory techniques.

All our products are made in France.

What is quantitative ATP-metry?

ATP-metry is a microbiological technique that evaluates the **overall bacterial load** of a water sample in less than **2 minutes**.



It is based on the detection of ATP molecules (energy-carrying molecule) which are found only in living cells:

« Any trace of ATP proves the existence of living organisms. »

▲ Watch the video

Once extracted from bacteria, the ATP reacts with our reagents, in particular with the luciferase enzyme which comes from the firefly's tail. The amount of light emitted is **proportional to the quantity of microorganisms**. It is measured by a luminometer.



ATP extracted
from bacteria



Enzyme reagents
DENDRIDIAG®

Bioluminescence reaction



Light emission (photons)
measured by a luminometer


Each measurement becomes **quantitative** because of the addition of a standard which takes into account environmental factors (temperature, pH, inhibitors...).

1 pg of ATP ≈ 1 000 bacteria

This easy and reliable on-site analysis is a great **decision support tool**. It allows you to validate treatments, to look for critical points on your facilities, manage crisis situations...

Applications


With GL BIOCONTROL's ATP-metry, you can:

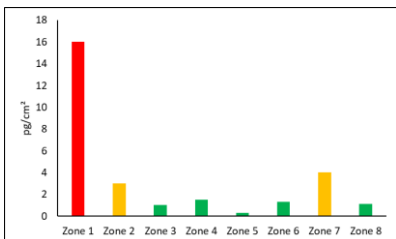

Control the facility cleanliness:

- Manage biofouling.
- Improve health risk management.


Avoid production shutdown.

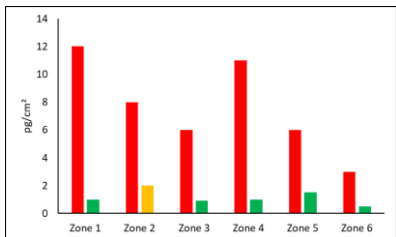

Assess operating procedure efficiency:

- Validate efficiency of:
 - Cleaning,
 - Disinfection.


Validate treatment strategy.

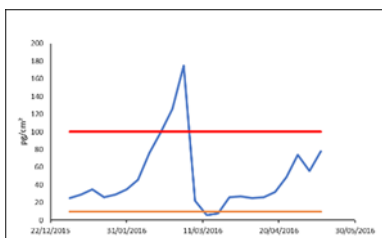

Manage biofilm formation in a pipe:

- Control and assess the fouling state of the pipe network.

Start and validate cleaning and disinfection procedures.

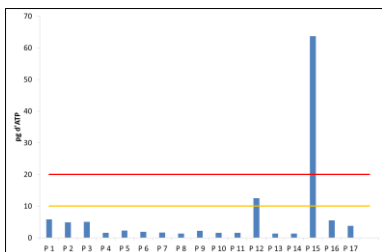

Identify the critical areas:

- Determine critical areas with important biomass growth.
- Adapt cleaning and disinfection strategy.

Highlight malfunctions.

Equipment needed

Reagent kit (50 tests)*

Product	Quantity
Dropper bottle EXTRACTANT	2
Dropper bottle STANDARD	1
Aluminium bag of 10 test tubes LUMITUBE	5

Consumable kit (50 tests)

Product	Quantity
Sterile swabs	50

Equipment

Product	Quantity
Stainless steel sampling template 20 cm ²	1
Luminometer KIKKOMAN SMART or equivalent	1
Tube holder	1

** The reagents can be stored at room temperature and in the dark for 3 months. To ensure an optimum conservation over 1 year, we advise you to keep them refrigerated (3 - 8°C).*

Protocole

Installation



Take the dropper bottles of **EXTRACTANT** and **STANDARD** out as well as a **LUMITUBE** pouch.



Make sure they are at room temperature (> 18°C) before use.

Take out the swabs as well as the sampling template. Turn on the luminometer and wait 8 seconds for the device calibration.

Sampling



Open a bag of **LUMITUBE** and take one lumitube.

Remove the cap with the tube holder, and place the lumitube on the rack supplied.



Get the swab out of its packaging being careful not to touch the cotton tip with your fingers or with the lab bench.


Squeeze 2 drops of **EXTRACTANT** on the tip and put the swab back in its packaging.



Clean the sampling template and place it on the surface.

Thoroughly swab the 20cm² surface. Pass at least two times on the surface on two different directions.

Measurement

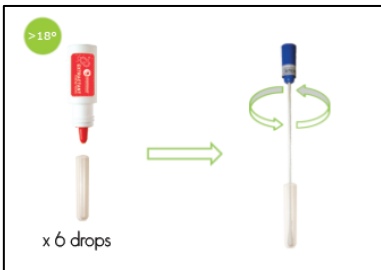

Squeeze 6 drops of **EXTRACTANT** in the lumitube.

Insert the swab in the lumitube and homogenize correctly by turning the swab using the rod.



Break off the top of the swab while maintaining it in the lumitube.


Fix the tube to its holder. Place them in the luminometer and press ENTER to start the measurement.

Write down the R1 result (in RLU).



Immediately, get the tube holder with the lumitube out of the luminometer.

Take the lumitube off the tube holder and add one drop of **STANDARD** in the middle of the lumitube.

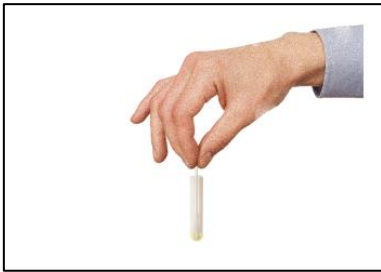

Correctly homogenize the mix by turning the swab inside the test tube to get a good calibration.



Fix the tube to its holder.

Place them in the luminometer and press ENTER to start the measurement.

Write down the R2 result (in RLU).

Interpretation of results

To interpret the results, you have two choices:

- The **DENDRIDIAG® APP** for smartphone and tablet,
- The Excel file.



Enter the surface sampled, R1 and R2 values measured by the luminometer.

Calculations are made automatically and are given in:

- **picogram** per square centimeter (pg ATP/cm²),
- total flora in **equivalent bacteria** per square centimeter (eq.bact./cm²)
- total flora in **logarithm** per square centimeter (LOG eq.bact./cm²)

Thanks to a color code, you are informed if a corrective action is required.

The Excel file provides traceability of results and allows you to monitor the level of bacteria across time and space.

[illegible]

If you **monitor your facility periodically**, fill in the « BIOMONITORING » sheet. A graph is automatically generated in the « GRAPH BIOMONITORING » sheet.


⇒ *It is advisable to monitor, at least once a week, the critical areas.*

If you **do a mapping of your facility**, fill in the « CARTOGRAPHY » sheet. A graph is automatically generated in the « GRAPH CARTOGRAPHY » sheet.


⇒ It is advisable to perform a measurement before and after cleaning or disinfection procedure for each critical area.

Warning and alarm thresholds were established based on our experience. These thresholds should be refined based on the first results obtained on your facility. The following arrows will help you **interpret the results**:


Operating swimming pools (in LOG eq.bact./cm²):

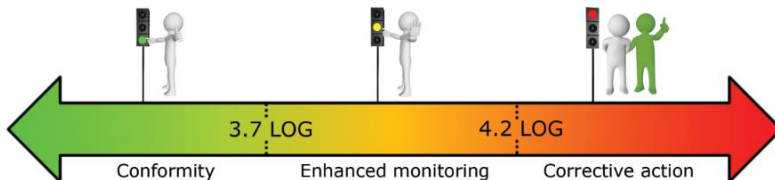

Surfaces of food processing industry (in pgATP/cm²):

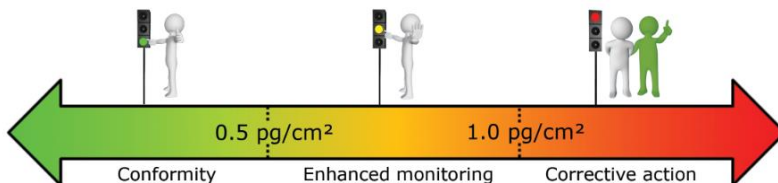

Surfaces of water pipes (in pgATP/cm²):

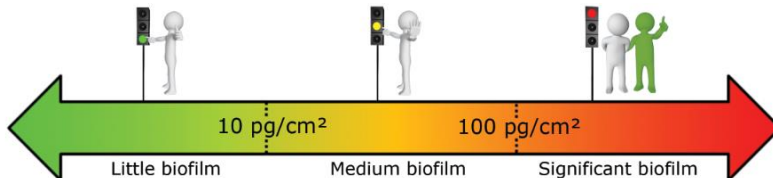

We consider that when a measurement is:

- **Below the warning threshold**, the facility is under control,
- **Between the warning and the alarm threshold**, the facility does not present an immediate biohazard. A corrective action is recommended if 3 consecutive measurements are above the warning threshold,
- **Above the alarm threshold**, the facility is not under control. A quick corrective action is recommended.

In the Excel file, the result is displayed in green, orange or red depending if it is under the warning threshold, between the warning and the alarm threshold or above the alarm threshold.

Troubleshooting

Problem	Possible cause and correction
"Low sensitivity. If necessary, sample a larger surface." displayed in the Excel file.	<p>The enzyme contained in the lumitube is not sufficiently active (out-of-date or degraded) or the EXTRACTANT is too cold.</p> <p>Warm up the reagent to a temperature above 18°C and filter a larger volume of water. If the problem remains, perform a Control of the reagent efficiency (cf. page 14).</p>
"Control the mixing of the STANDARD 1000, the temperature and condition of the reagents." displayed in the Excel file.	<p>Standardization was not successful. Tap the bottom of the lumitube on a flat surface, homogenize and restart the measurement.</p> <p>If the problem remains, the enzyme contained in the lumitube is not sufficiently active (out-of-date or degraded) or the EXTRACTANT is too cold. Warm up the reagent and perform Control of the reagent efficiency (cf. page 14).</p>
"Sample highly contaminated. If necessary, sample a smaller surface." displayed in the Excel file.	<p>The ATP concentration of the sample is too high. Restart the analysis by sampling a smaller surface.</p>

Controls



To ensure proper functioning of your luminometer, we advise you to conduct an annual maintenance. We remain at your disposal to assist you.

Control of the luminometer contamination

- Place the tube holder in the luminometer and press ENTER,
- The result should be less or equal to 2 RLU. If not, with a cotton swab, wipe the internal surfaces of the measurement chamber.

Control of the reagent contamination

- In a lumitube, put 4 drops of **EXTRACTANT**,
- Fix the lumitube to its holder,
- Place it in the luminometer and press ENTER,
- The result should be less or equal to 5 RLU. If not, your dropper bottle of **EXTRACTANT** is probably contaminated. Contact GL Biocontrol for technical support.

Control of the reagent efficiency

- In a lumitube, put 4 drops of **EXTRACTANT** and 1 drop of **STANDARD** (reagent temperature must be above 18°C),
- Fix the lumitube to its holder,
- Properly homogenize the lumitube,
- Place it in the luminometer and press ENTER,
- For a good efficacy of the reagents, the result should be higher than 400 RLU. If not, the lumitubes are probably degraded. Discard the bag of **LUMITUBE** and choose a new one.

Contact

For further information or assistance on result interpretation, on the protocol or for commercial information, contact by email or by phone:

Yannick FOURNIER

Sales engineer

5, avenue de l'Europe - Hélioparc

34 830 CLAPIERS (FRANCE)

Phone: +33 (0)9 67 39 35 20

Phone: +33 (0)6 33 64 42 29

Email: y.fournier@gl-biocontrol.com

Clément FAYE

Research engineer

5, avenue de l'Europe - Hélioparc

34 830 CLAPIERS (FRANCE)

Phone: +33 (0)9 67 39 35 20

Phone: +33 (0)6 72 70 46 98

Email: c.faye@gl-biocontrol.com

A tutorial video is available on the flash drive supplied with the luminometer or on our website (<https://www.gl-biocontrol.com/en/video-tutorial>), or here:



◀ Video tutorial

3 easy ways to order

@ by e-mail at sales@gl-biocontrol.com,

☎ by phone at + 33 (0) 9 67 39 35 20,

✉ by mail at GL BIOCONTROL - 5, avenue de l'Europe, Hélioparc - 34,830 CLAPIERS (FRANCE).

>>> www.gl-biocontrol.com

GLBIOCONTROL

R A P I D M I C R O B I A L D I A G N O S T I C

5, avenue de l'Europe – Hélioparc

34 830 CLAPIERS (FRANCE)

Phone: +33 (0)9 67 39 35 20

Email: contact@gl-biocontrol.com

Web: www.gl-biocontrol.com